



SEP 06, 2022

OPEN  ACCESS

**DOI:**  
[dx.doi.org/10.17504/protocols.io.3byl4je4rlo5/v1](https://dx.doi.org/10.17504/protocols.io.3byl4je4rlo5/v1)

**External link:**  
<https://www.biorxiv.org/content/10.1101/2022.02.18.481043v1.full>

**Protocol Citation:** Itika Saha, F. Ulrich Hartl, Mark S. Hipp 2022. Detection of Tau ubiquitylation. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.3byl4je4rlo5/v1>

**MANUSCRIPT CITATION:**  
 The AAA+ chaperone VCP disaggregates Tau fibrils and generates aggregate seeds  
 Itika Saha, Patricia Yuste-Checa, Miguel Da Silva Padilha, Qiang Guo, Roman Körner, Hauke Holthusen, Victoria A. Trinkaus, Irina Dudanova, Rubén Fernández-Busnadiego, Wolfgang Baumeister, David W. Sanders, Saurabh Gautam, Marc I. Diamond, F. Ulrich Hartl, Mark S. Hipp  
 bioRxiv 2022.02.18.481043;  
 doi:  
<https://doi.org/10.1101/2022.02.18.481043>

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
## Detection of Tau ubiquitylation

Itika Saha<sup>1</sup>, F. Ulrich Hartl<sup>2,3</sup>, Mark S. Hipp<sup>2,3</sup>

<sup>1</sup>Department of Cellular Biochemistry, Max Planck Institute of Biochemistry, Am Klopferspitz 18, 82152 Martinsried, Germany;

<sup>2</sup>Department of Biomedical Sciences of Cells and Systems, University Medical Center Groningen, University of Groningen, Antonius Deusinglaan, 1, 9713 AV Groningen, The Netherlands;

<sup>3</sup>School of Medicine and Health Sciences, Carl von Ossietzky University Oldenburg, 26129 Oldenburg, Germany



Felix Kraus

### ABSTRACT

This protocol describes the detection of ubiquitylated Tau from HEK293 cells stably expressing and propagating aggregates of Tau repeat domain fused to YFP (Sanders et al. Neuron, 2014; Saha et al, BioRxiv, 2022).

**License:** This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

**Protocol status:** Working  
We use this protocol and it's working

**Created:** Sep 05, 2022

**Last Modified:** Dec 26, 2022

**PROTOCOL integer ID:**  
69589

**Keywords:** ASAPCRN

- 1 Harvest cells from 2 confluent wells of a 6 well plate.
- 2 Lyse aggregate-containing cells by vortexing in cold RIPA buffer (Thermo) supplemented with protease inhibitor cocktail (Roche) and DNase. 20 mM N-ethylmaleimide should be included in the lysis buffer to inhibit the activity of deubiquitinating enzymes.
- 3 Briefly sonicate lysate and centrifuge at 2,000 x g for 5 min to remove cell debris. 5m
- 4 Collect supernatant, determine protein concentration and normalize across samples.
- 5 Dilute 1 mg protein in a total volume of 600 µL RIPA buffer.

- 6 Add 50  $\mu$ L anti-GFP bead slurry ( $\mu$ MACS GFP Isolation kit, Miltenyi Biotec) to diluted lysate.
- 7 Incubate for 1 h at 4 °C in a rotating wheel at 10 rpm. 1 h
- 8 Before the end of 1 h, place  $\mu$ -columns (Miltenyi Biotec) in the magnetic field of a  $\mu$ MACS Separator (Miltenyi Biotec) and equilibrate columns by applying 250  $\mu$ L RIPA buffer. Allow complete flow-through.
- 9 Apply cell lysates and beads to  $\mu$ -columns. Allow complete flow-through.
- 10 Wash columns 4 times with 1 mL 0.1% SDS/PBS.
- 11 Elute by applying 50  $\mu$ L pre-heated (95 °C) 1x SDS sample buffer.
- 12 Analyze eluates by immunoblotting with antibodies against GFP or ubiquitin.