



SEP 06, 2022

OPEN ACCESS

DOI:
dx.doi.org/10.17504/protocols.io.x54v9den4g3e/v1

External link:
<https://www.biorxiv.org/content/10.1101/2022.02.18.481043v1.full>

Protocol Citation: Miguel Da Silva Padilha, Irina Dudanova, Itika Saha, F. Ulrich Hartl, Mark S. Hipp 2022. Immunohistochemistry of rTg4510 mouse brain. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.x54v9den4g3e/v1>

MANUSCRIPT CITATION: Itika Saha, Patricia Yuste-Checa, Miguel Da Silva Padilha, Qiang Guo, Roman Körner, Hauke Holthusen, Victoria A. Trinkaus, Irina Dudanova, Rubén Fernández-Busnadiego, Wolfgang Baumeister, David W. Sanders, Saurabh Gautam, Marc I. Diamond, F. Ulrich Hartl, Mark S. Hipp bioRxiv 2022.02.18.481043; doi: <https://doi.org/10.1101/2022.02.18.481043>

🌐 Immunohistochemistry of rTg4510 mouse brain


Miguel Da Silva Padilha¹, Irina Dudanova¹, Itika Saha², F. Ulrich Hartl², Mark S. Hipp^{3,4}

¹Molecular Neurodegeneration Group, Max Planck Institute of Neurobiology, 82152 Martinsried, Germany;

²Department of Cellular Biochemistry, Max Planck Institute of Biochemistry, Am Klopferspitz 18, 82152 Martinsried, Germany;

³Department of Biomedical Sciences of Cells and Systems, University Medical Center Groningen, University of Groningen, Antonius Deusinglaan, 1, 9713 AV Groningen, The Netherlands;

⁴School of Medicine and Health Sciences, Carl von Ossietzky University Oldenburg, 26129 Oldenburg, Germany

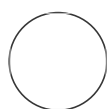

Felix Kraus

ABSTRACT

This protocol describes immunohistochemical staining of the Tau transgenic mice rTg4510 (Santacruz et al, 2005) brain section for Tau aggregates phosphorylated at Ser202/Thr205 and the AAA+ ATPase Valosin-containing protein (VCP). Experiments involving animal models must be performed in accordance with relevant institutional guidelines and regulations.

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working
We use this protocol and it's working

Created: Sep 05, 2022

Last Modified: Dec 26, 2022

PROTOCOL integer ID:
69594

Keywords: ASAPCRN

1 Deeply anesthetize mice with 1.6% Ketamine/0.08% Xylazine and transcardially perfuse with PBS followed by 4% paraformaldehyde (PFA)(Santa Cruz) in PBS.

2 Dissect brains out of the skull and post-fix in 4% PFA in PBS overnight.

12h

3 Embed fixed tissue in agarose and section into 40 µm thick sections using a vibratome (VT1000S, Leica).

Note

NOTE: The sections can be stored in PBS 0.05% sodium azide at 4°C.

4 Permeabilize sections with 0.5% Triton X-100 and wash with PBS.

5 Incubate sections in blocking solution consisting of 0.2% BSA (w/v), 5% donkey serum (v/v)

30m

(abcam), 0.2% lysine (w/v) (Sigma-Aldrich), 0.2% glycine (w/v) (Sigma-Aldrich) in PBS for 30 min at room temperature (RT).

- 6** Incubate sections with primary antibodies (anti-phospho-Tau AT-8 (Thermo, Cat# MN1020, 1:300); anti-VCP (Novus Biologicals, Cat# NB 100-1558, 1:500) diluted in 0.3% Triton X-100 (v/v), 2% BSA (w/v) in PBS overnight at 4°C. 12h
- 7** Wash sections in PBS and incubate with secondary antibodies and Neurotrace 640/660 (ThermoFisher, Cat# N21483, 1:500) diluted in 0.3% Triton X-100, 3% donkey serum (v/v) for 2 h at RT. 2h
- 8** Stain nuclei with 0.5 µg/ml DAPI.
- 9** Mount sections on Menzer glass slides using Prolong Glass fluorescence (Invitrogen) mounting medium.